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울산과학기술원 연구지원본부 기기분석실  
UNIST Central Research Facilities  
UNIST Materials Characterization Lab

## CONTACT

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# SEM Data 이동

The screenshot shows a Windows 10 desktop with a blue background. On the left, there are three numbered steps: 1) Open Chrome, Internet Explorer, or Firefox (Windows XP). 2) Type 10.24.9.32 (Online, Windows 10) or 100.100.100.30 (Offline, Windows 10 <). On the right, there is a section titled '3) Login UCRFSERVER' with a login form. The form has a title 'UCRFSERVER', a user ID field, a password field, a '로그인 유지' checkbox, and a '로그인' button. Below the instructions, there are two browser window screenshots. The top one is labeled 'Online' and shows the address bar with '10.24.9.32'. The bottom one is labeled 'Offline' and shows the address bar with '100.100.100.30'. The Windows taskbar is visible at the bottom.

1) Open  
Chrome  
Internet Explorer  
Firefox (Window XP)

2) Type  
10.24.9.32 (Online, Windows 10)  
100.100.100.30  
(Offline, Windows 10 <)

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Daily checklist in lab.

Lab title	Electron Microscopy Preparation	Bldg./NO.	102-B115	Date	
Division	checkpoint			check	
General Safety	Laboratory overall cleanliness condition.				
	Smoking or bringing food into the lab.				
	Management status of experimental equipment such as safety regulations, safety signs, personal protective equipment, first aid, etc.				
	Checking the presence of a pre-hazard risk analysis report.				
Electric	Checking the power supply status of unused electrical equipment and checking for overloaded outlets.				
	Using grounded outlets, checking damage on the insulating coating of electric wiring, electric wiring arrangement.				
	Checking ground conditions for preventing external or static disturbances of the instrument.				
	Non-load status around electric panelboards.				
Fire	Fire extinguisher sign, proper fire extinguisher and regular inspection status.				
	Emergency exits, escape routes, and any obstacle blocking the passage.				
	Storage of foreign substances around fire hydrant and fire extinguishers.				
Gas	Outdoor storage of gas containers, no risk of tipping over, and checking ventilation conditions.				
	Corrosion, deformation, nozzle lock status on the exterior of the gas containers and checking the packing time limit of gas containers.				
	Checking installation and operational status of gas leakage detection alarm, anti-backflow/anti-backfire prevention devices, neutralizing decontamination devices.				
	Attachment of pipe marks, gas facility boundaries/warning marks and operation status of regulators and valves.				
	Safe separation distance from surrounding fire hazards.				
Chemical	Keeping hazardous factors handling and management registers and MSDS.				
	Categorizing chemicals by description and storing chemical reagents in safety cabinets.				
Equipment check	SU8220 Cold FE-SEM user: Check the pressure of nitrogen gas				
Confirm	Inspector Signature				(sign)
	Lab Director Signature				(sign)

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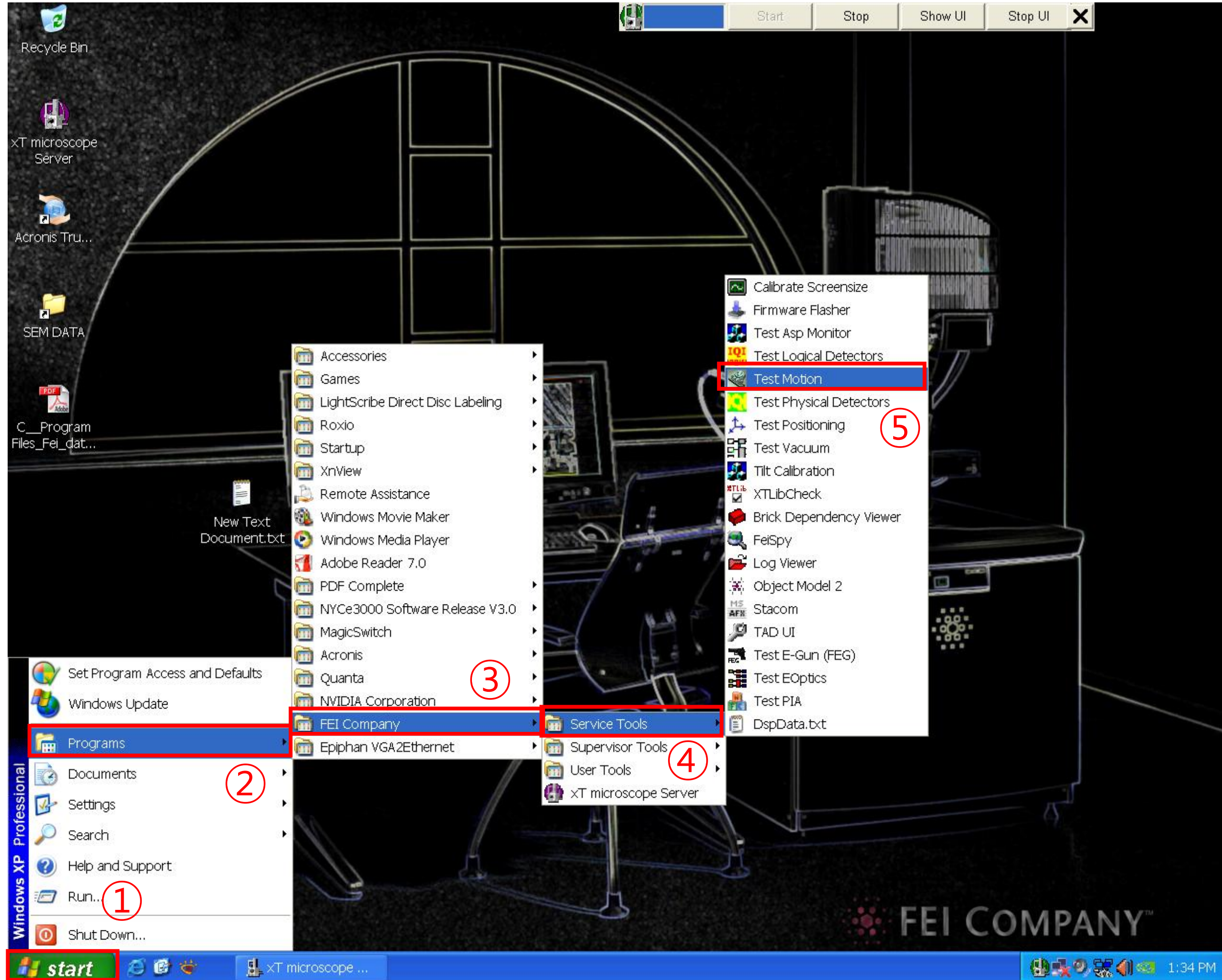
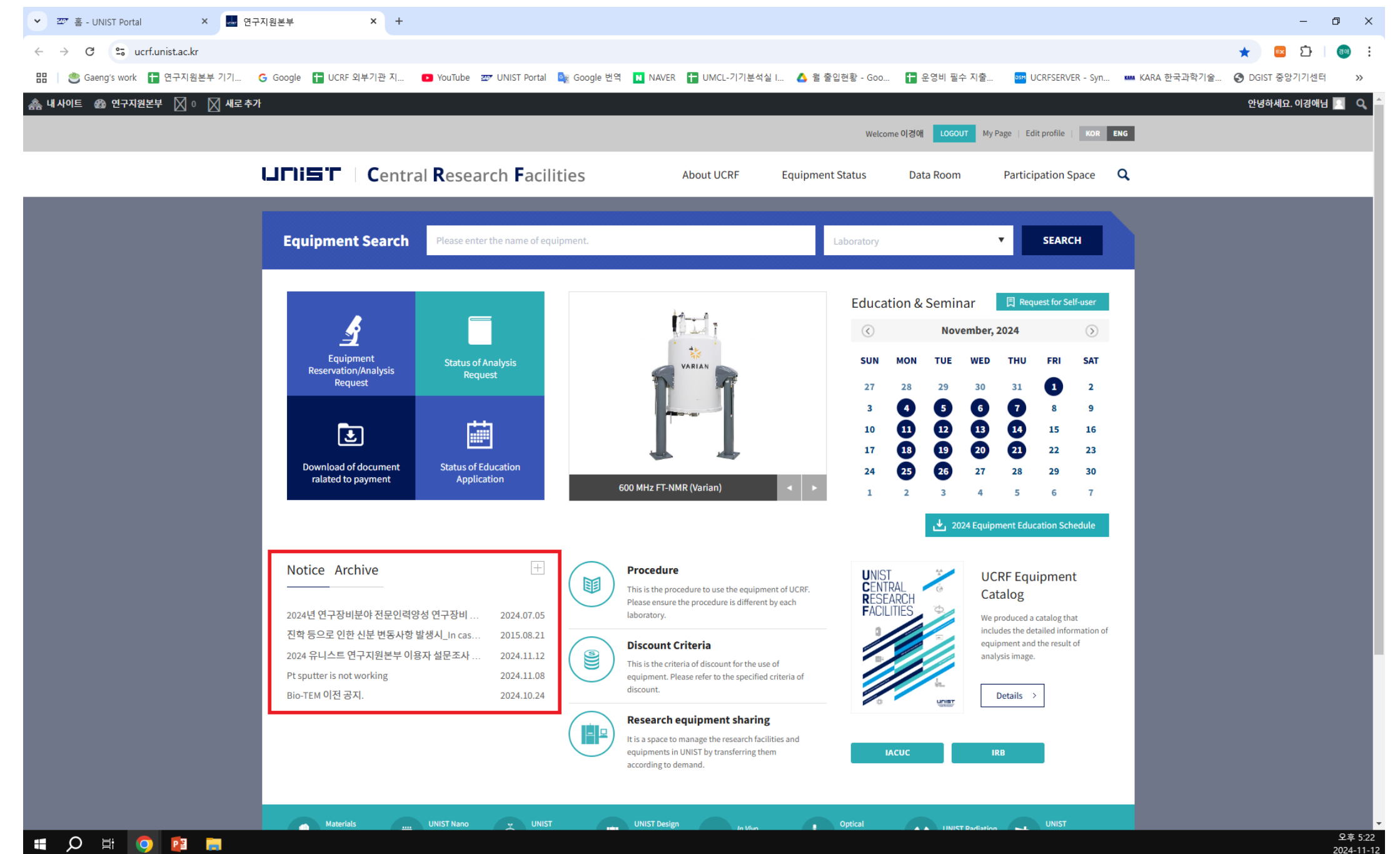
# Energy Dispersive X-ray Spectroscopy

The screenshot displays the EDAX software interface. At the top left is the EDAX logo with the tagline 'Smart Insight' and the AMETEK logo with 'MATERIALS ANALYSIS DIVISION'. The website address 'www.edax.com' is centered at the top. Below this is a 'Crystal Structures' panel with icons for Cubic (face centered, body centered), Hexagonal, Monoclinic, Orthorhombic, Tetragonal, and Rhombohedral. The main part of the interface is a periodic table where elements are color-coded by groups: IA (yellow), IIA (purple), IIIA (orange), IVA (green), VA (red), VIA (blue), VIIA (light blue), and VIIIA (dark blue). A legend at the bottom left identifies categories like Alkali Earth, Alkaline Earth, Rare Earth, Other Metals, Non-Metals, Halogens, Transition Metals, Metalloids, Noble Gases, and Actinides. The periodic table includes atomic numbers, symbols, and names for all elements from Hydrogen to Oganesson.

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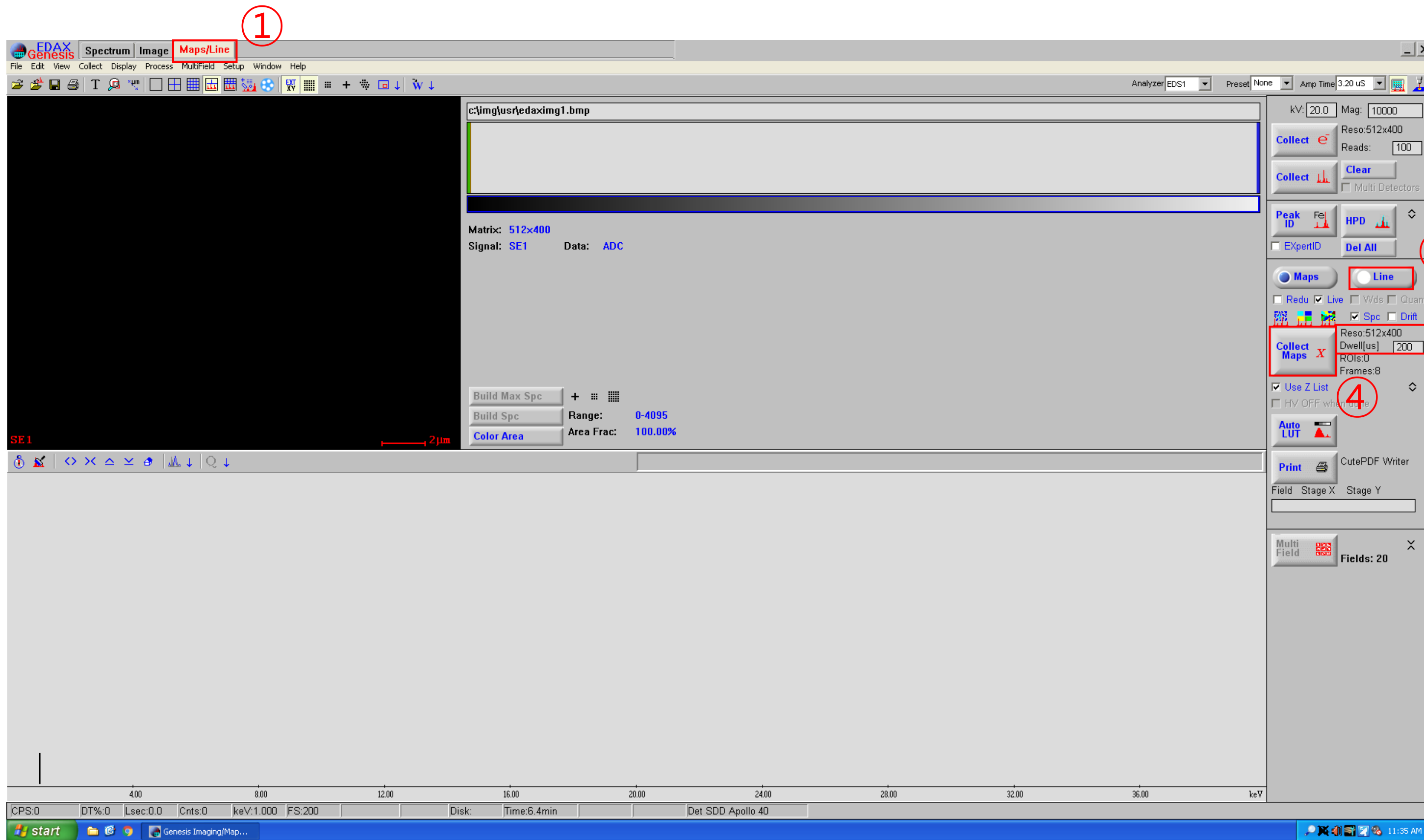
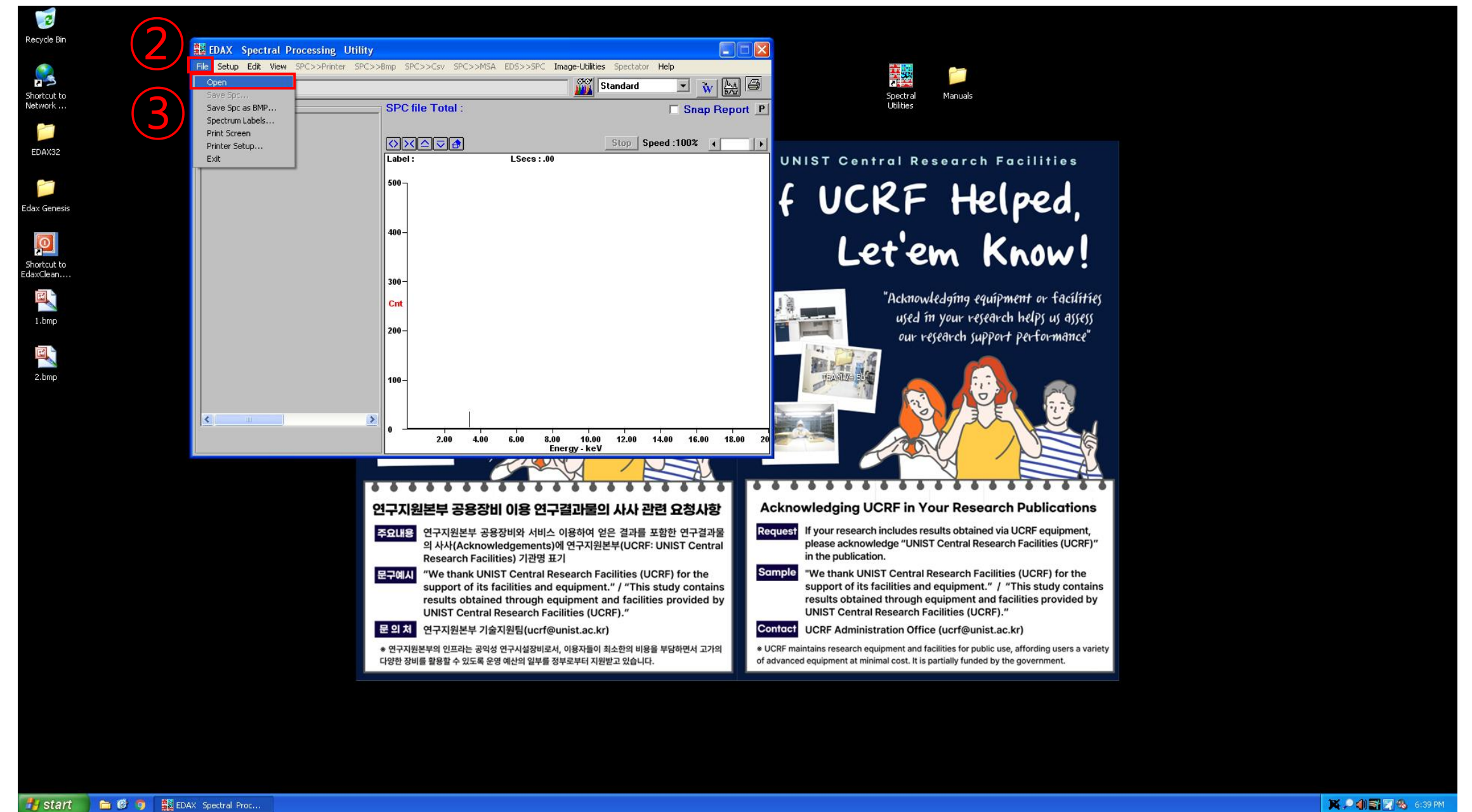
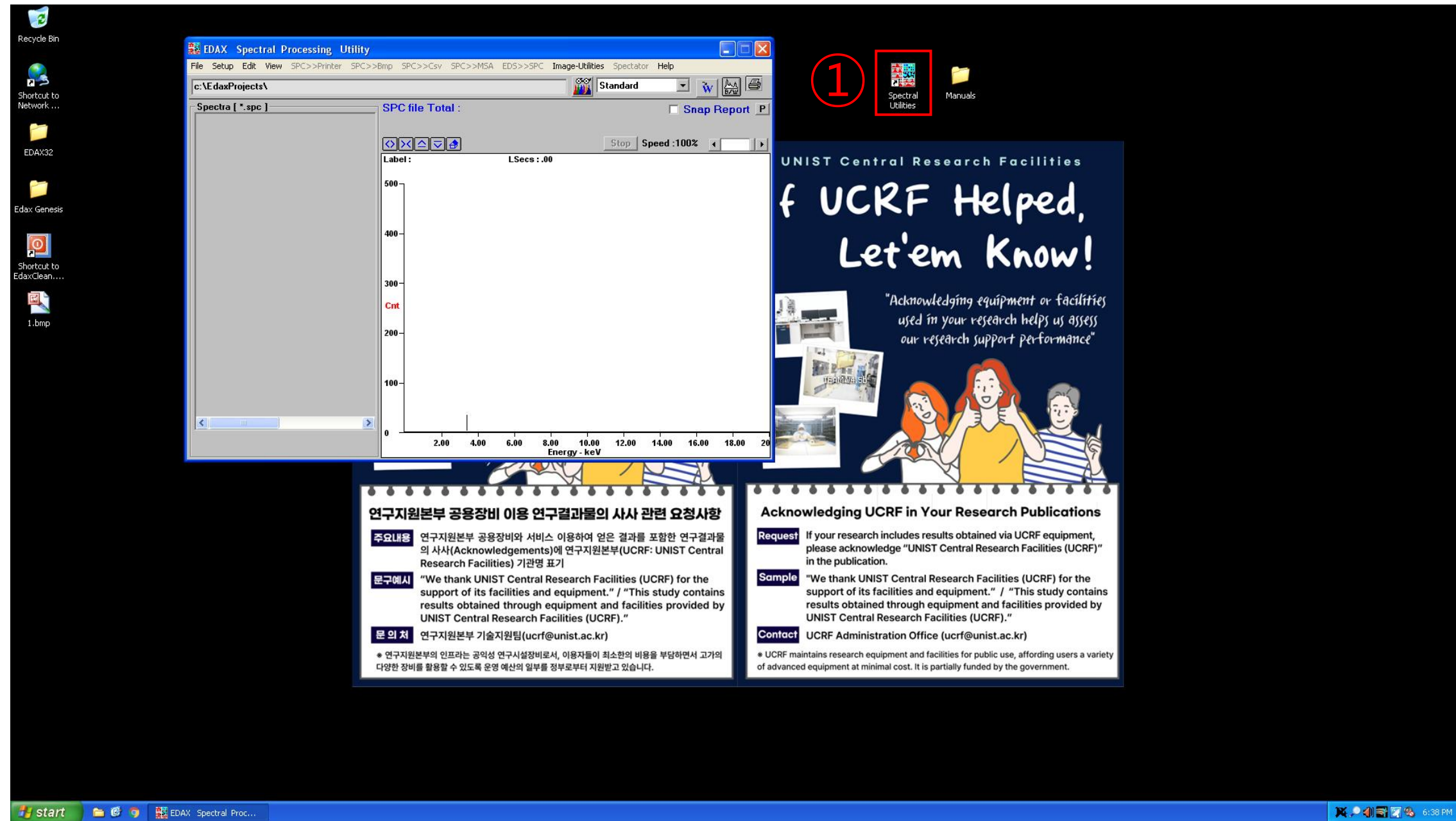
※ If the problem cannot be resolved, please contact the person in charge.

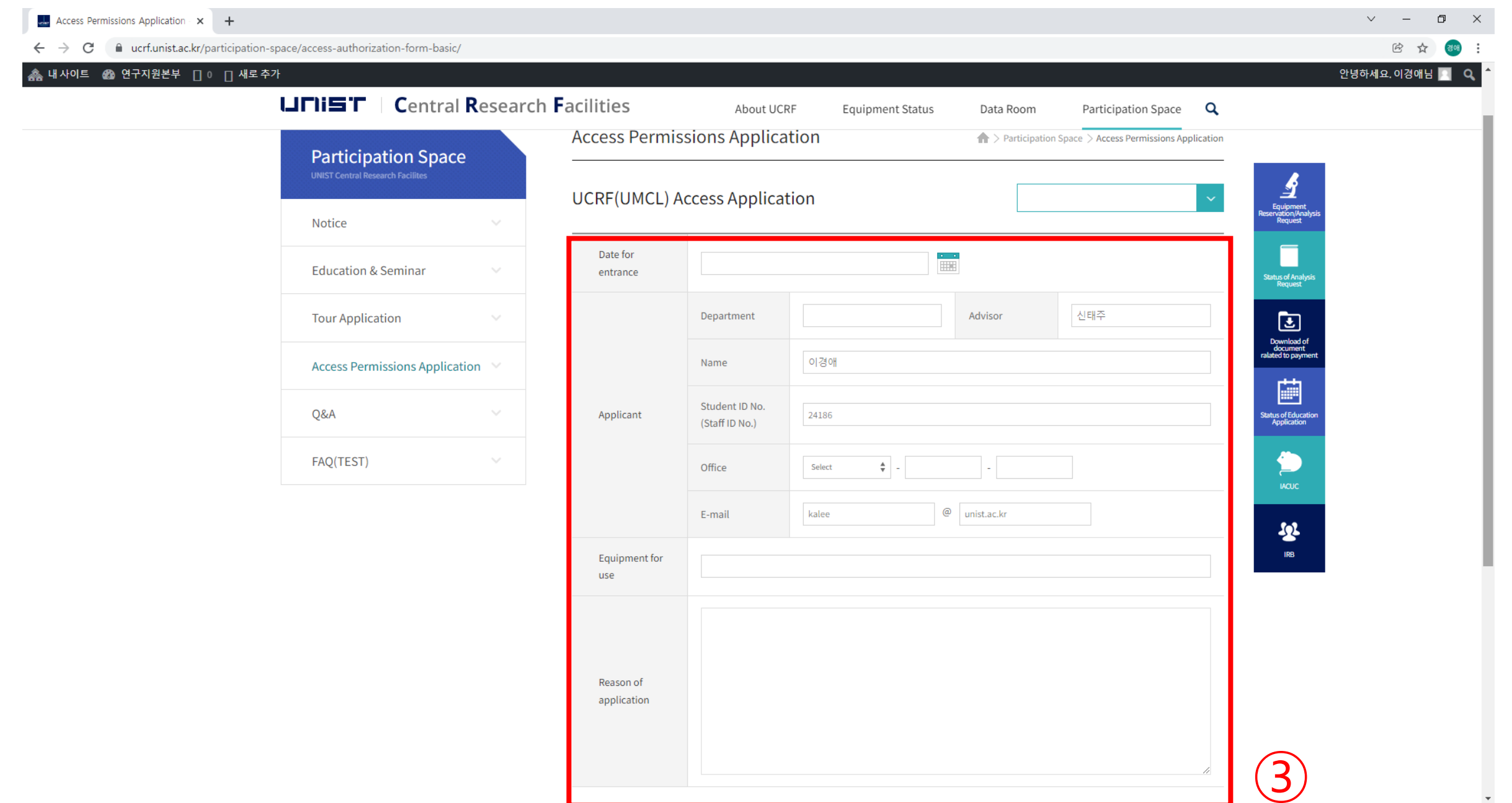
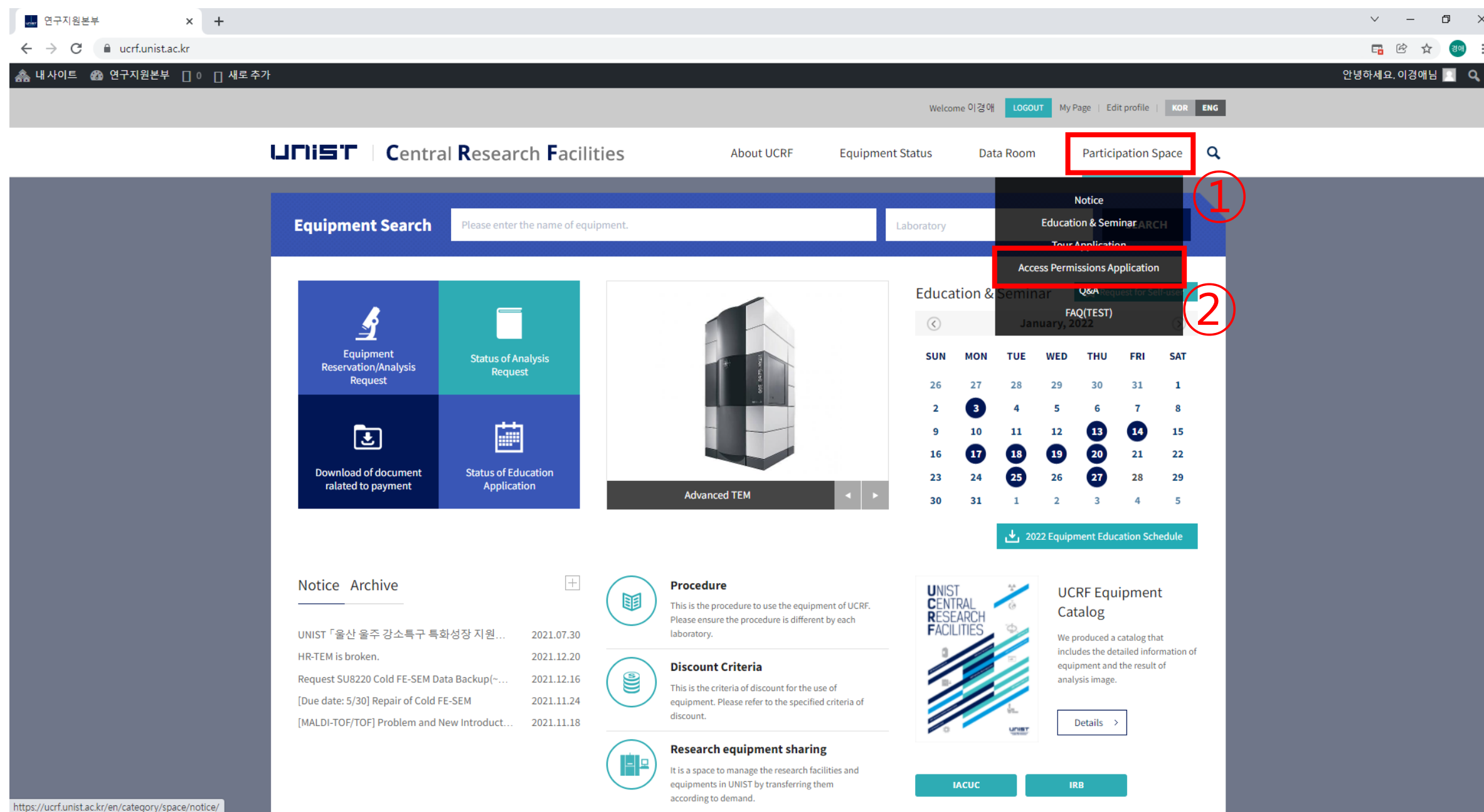
## Subject: Image Collection problem

Problem: No image collected (Black, Gray or White) from the microscope in Genesis. Image collection is okay in the EDAM or SG Shell.

If the image collection in the SG or EDAM Shell is okay, but the image collection in TEAM or Genesis is not okay, it may be due to a corrupted configuration file.

- ✓ Close all EDAX software
- ✓ Go to the **My Documents** folder and delete all **gen\*.ini** files such as **geneds.ini**, **genimg.ini**, **genpart.ini**, etc.
- ✓ Check if image collection in Genesis works.
  
- ✓ If the above does not work do the next steps:
- ✓ Close all EDAX software
- ✓ Use Windows Explorer to browse the c:\ drive
- ✓ Go to c:\edax32\sys folder
- ✓ Locate and then double click on **edaxclean.exe**
- ✓ Go to c:\edax32\img folder
- ✓ Delete file **edi32coll.cfg**.
- ✓ Delete the file **gpuscan.cfg**
- ✓ Reboot the pc.
- ✓ Check the image collection in Genesis
  
- ✓ If the above does not work do the next steps:
- ✓ Close all EDAX software
- ✓ Use Windows Explorer to browse the c:\ drive
- ✓ Go to c:\edax32\Genesis folder
- ✓ Select and **RENAME: Genesis.ini** to **Genesis.old**
- ✓ Select and **RENAME: Preference.ini** to **Preference.old**
- ✓ Check the image collection in Genesis

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